

| CSOP Title | |
|------------------------|--|
| CSOP No. | CSOP60 |
| Version | 1.1 |
| Location | UQ Node/Centre-wide |
| Policy/Procedure Link | UQ- Equipment |
| Risk Assessments | |
| Approved by | <i>Lindsay Shaw</i> |
| Date Approved | 13/10/2021 |
| Date Effective | 05/08/2021 |
| Next Review Date | 05/08/2022 |
| Contact for Assistance | plantsuccess@uq.edu.au |

1.0 Scope

To outline the use of the growth cabinets. This includes cleaning, what to do if it malfunctions/breaks down.

2.0 Prescribed Actions

- Open door
- Put plants into the cabinet
- Close the door (and watch fingers)
- Make sure you clean up after each experiment
 - Remove seeds with brush or vacuum
 - Clean with ethanol or F10
- To change conditions, please ask a postdoc for help as the program is complicated to use.
- If the cabinet breaks down, contact number on the cabinet.
- If any disease or insects are noticed on any plants, notify other users of the cabinet and decide on best course of action to eradicate. Do not use

LABORATORY STANDARD OPERATING PROCEDURE (LSOP)

ARC COE for Plant Success in Nature and Agriculture: *Using Growth Cabinets*

other cabinets and growth rooms after opening contaminated cabinet to prevent spread of disease.